




BATTERY HOLDERS FOR SWAGELOK

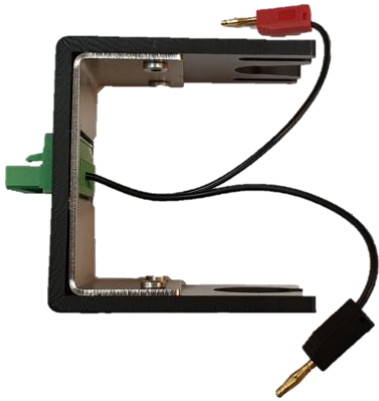

2 Electrodes

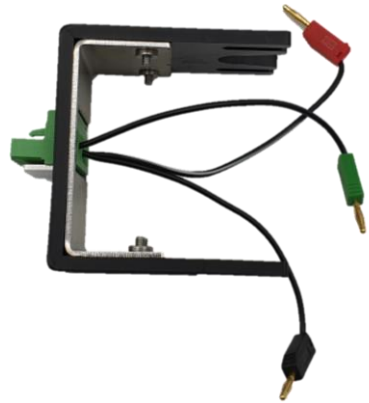


3 Electrodes

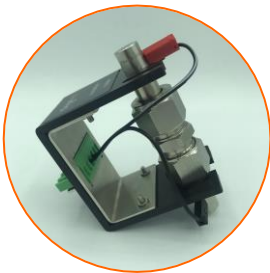

Battery Holder

HOW IT WORKS

1 Insertion of the Swagelok

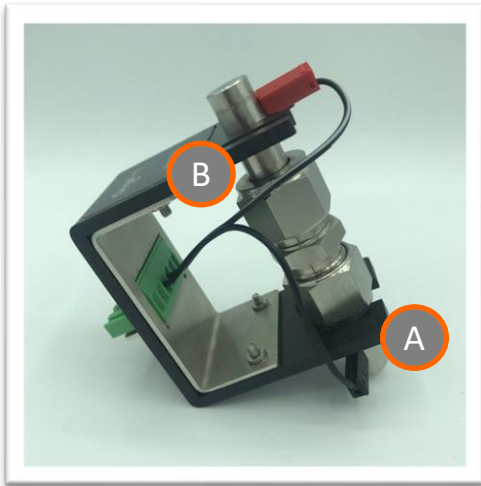


Insert the Swagelok in the Battery Holder from the bottom part (A). Then, "clips" the rod in the top part (B).

TIPS: To be sure the Swagelok is well inserted; you can check its stability. If well inserted into the Battery Holder, the Swagelok must not move.

Once inserted, connect as below:


- Red connector (**WRK**) to the Anode.
- Black connector (**AUX**) to the Cathode.

2 Connection to the potentiostat

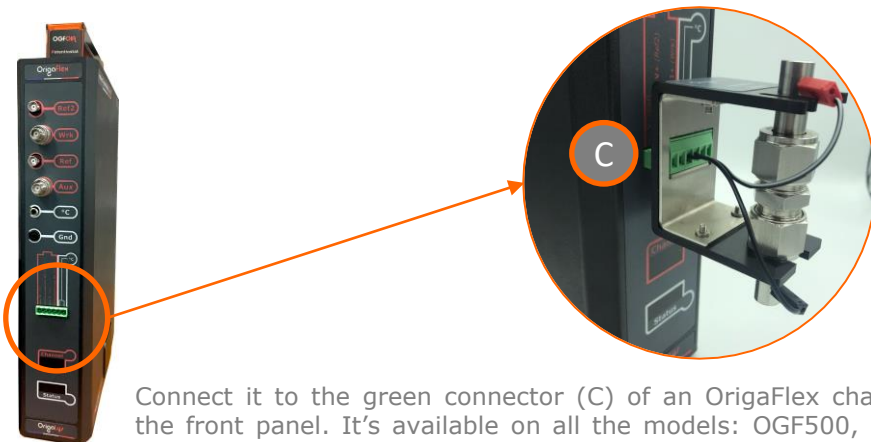


Connect it to the green connector (C) of an OrigaFlex channel, on the front panel. It's available on all the models: OGF500, OGF01A, OGF05A and OGF10A.

Swagelok 2 Electrode Type

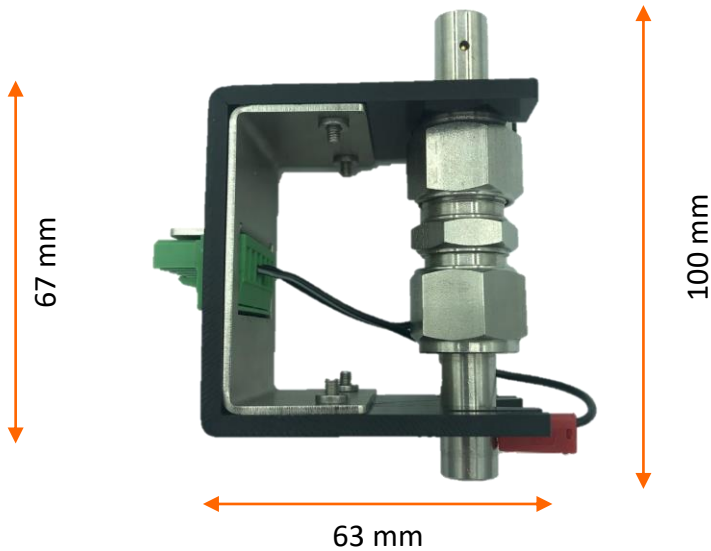

90 g

Only the Battery Holder

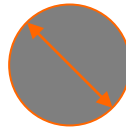

12 mm

Internal diameter

Other Battery Holders



Swagelok

✓ AR00884



For long cable (3m)

✓ CAD051RV1




For coin cell CR2032

✓ AR01190

Battery Holder

HOW IT WORKS

1 Insertion of the Swagelok

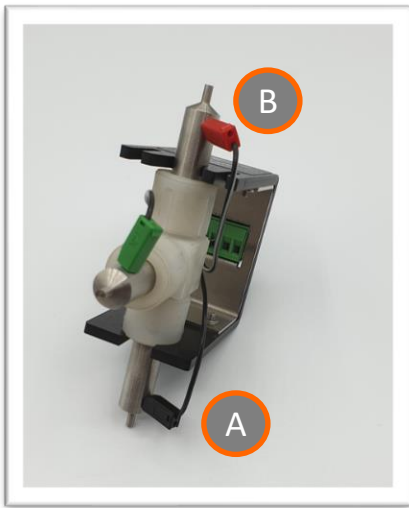


Insert the Swagelok in the Battery Holder from the button part (A). Then, "clips" the rod in the top part (B).

TIPS: To be sure the Swagelok is well inserted; you can check its stability. If well inserted into the Battery Holder, the Swagelok must not move.

Once inserted, connect as below:


- Red connector (**WRK**) to the Anode.
- Black connector (**AUX**) to the Cathode.
- Green connector (**REF**) for the Reference.

2 Connection to the potentiostat

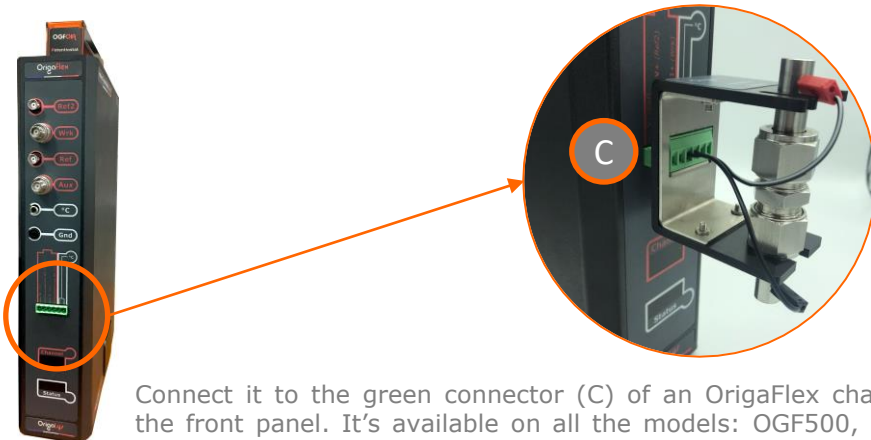


Connect it to the green connector (C) of an OrigaFlex channel, on the front panel. It's available on all the models: OGF500, OGF01A, OGF05A and OGF10A.

Swagelok 3 Electrode Type

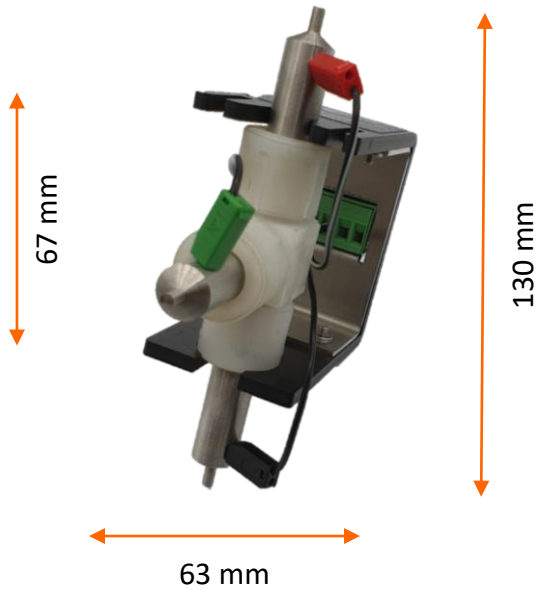

158 g

Only the Battery Holder

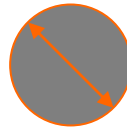

12 mm

Internal diameter

Other Battery Holders



Swagelok

✓ AR00885



For long cable (3m)

✓ CAD051RV1



For coin cell CR2032

✓ AR01190

BATTERIES



www.origalys.com


CORROSION



www.origalys.com




ORIGAFLEX

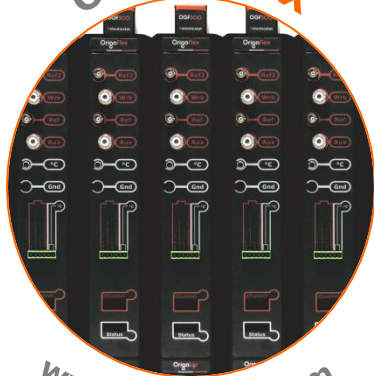

www.origalys.com

RESEARCH

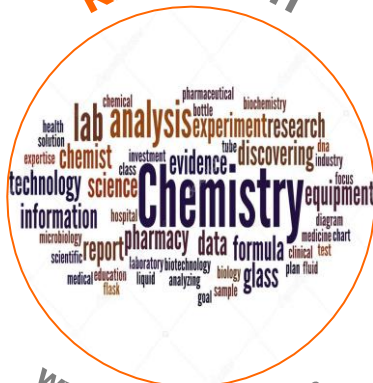

www.origalys.com

Subject to change
16-07-2021

Country Representative

OrigalyS ElectroChem SAS

Les Verchères 2
62A, avenue de l'Europe
69140 RILLIEUX-la-PAPE
FRANCE

+33 (0)9 54 17 56 03

+33 (0)9 59 17 56 03

contact@origalys.com